

78919969 imaging chamber – 5 mouse with Equaflow 5 manifold Quick-Start Guide

Items included:

Imaging chamber
Steel plate
Tubing connections

Equaflow 5 mouse manifold
Hepa filters (3)

- Remove all items from the packing material and inspect for any shipping damage. Notify Patterson immediately if any damage is noted.
- Remove the Equaflow 5 manifold from the inside of the chamber by turning the quick connects counter-clockwise. Install the dividers into the manifold as shown below and re-install back into the chamber:



- Connect the tubing pieces and Hepa filters as shown below (note the connections that will need to be made to the anesthesia system):

